


Installation instructions

1. Download APMALDI Firmware update
2. Unzip and run install.bat
3. Accept server host key if needed (see image bellow)

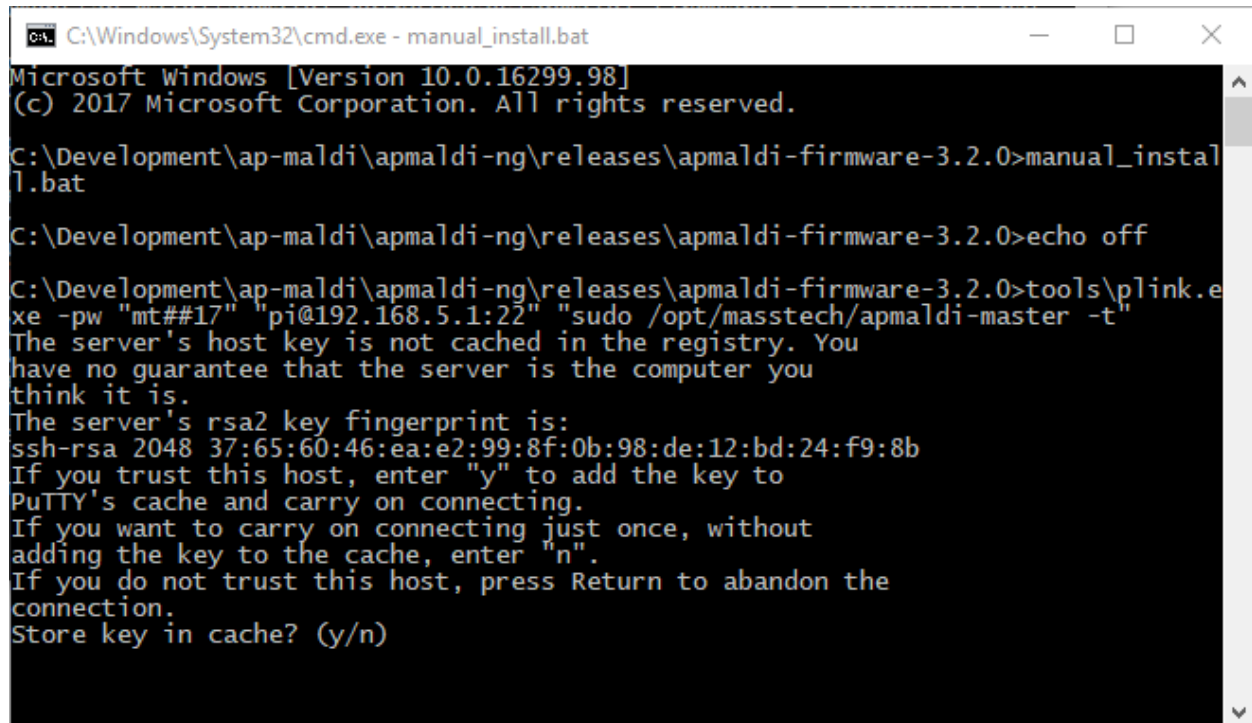

```
C:\Windows\System32\cmd.exe - manual_install.bat
Microsoft Windows [Version 10.0.16299.98]
(c) 2017 Microsoft Corporation. All rights reserved.

C:\Development\ap-maldi\apmaledi-ng\releases\apmaledi-firmware-3.2.0>manual_install.bat

C:\Development\ap-maldi\apmaledi-ng\releases\apmaledi-firmware-3.2.0>echo off

C:\Development\ap-maldi\apmaledi-ng\releases\apmaledi-firmware-3.2.0>tools\plink.exe -pw "mt##17" "pi@192.168.5.1:22" "sudo /opt/masstech/apmaledi-master -t"
The server's host key is not cached in the registry. You
have no guarantee that the server is the computer you
think it is.
The server's rsa2 key fingerprint is:
ssh-rsa 2048 37:65:60:46:ea:e2:99:8f:0b:98:de:12:bd:24:f9:8b
If you trust this host, enter "y" to add the key to
PuTTY's cache and carry on connecting.
If you want to carry on connecting just once, without
adding the key to the cache, enter "n".
If you do not trust this host, press Return to abandon the
connection.
Store key in cache? (y/n)
```